

- Fix with 2% Paraformaldehyde (~ 30 minutes)
- Permeabilize in PBS/ 0.01M Glycine? 0.1% Triton-X
  - 1 hour @ RT
- Wash in PBS
  - 3X, ~ 5 minutes each
- Block in 5% BSA/PBS
  - 1 hour @ RT
- Wash in PBS
  - 3X, ~ 5 minutes each
- Apply Primary Ab (1:250 in PBS)
  - Overnight @4 C
- Wash in PBS
  - 3X, ~ 5 minutes each

\*\*\*\*\***COVER WITH FOIL**\*\*\*\*\*

---

- Apply Secondary Ab (1:200 in PBS)
  - 1 hour @ 37 C
- Wash in PBS
  - 3X, ~ 5 minutes each
- Apply DAPI (1:1000) and Phalloidin 488 (1:250)
  - 1 hour @ RT
- Wash in PBS
  - 3X, ~ 5 minutes each
- Either image in well-plate on EVOS or Mount on slide with DABCO